

Flexible Arm Electrode Holder

TPS Part No 130088

Thank you for purchasing a TPS Flexible Arm Electrode Holder. This accessory will simplify the use of your instrument, and help protect electrodes against accidental breakage.

Unpacking Information

You should have received:

1. Electrode Support Arm
2. Electrode Support Base
3. Electrode Holder Clamp

If any of these items are missing, please contact TPS immediately.

Assembly

1. Insert the Electrode Support Arm into the Base. To firmly secure the arm, press it into the base until it locks into place.

NOTE: Do not twist the Electrode Support Arm when inserting into the base.

2. Slip the Electrode Holder Clamp onto the plastic rod at the end of the arm.

Operation

1. Fit up to two 12mm diameter electrodes into the large clips and one 6mm electrode into the small clip.
2. Position the arm and electrode holder as required. Tighten or loosen the thumb screws on the electrode to adjust tension as required.

Maintenance

1. The electrode arm may be lubricated at the joints from time to time. Use only silicon based lubricants.
2. Wipe clean with a moist, soft cloth.

TPS Pty Ltd
4 Jamberoo Street,
Springwood, Brisbane,
Australia, 4127

Phone	: (07) 32 900 400	Fax	: (07) 3808 4871
International	: 61 7 32 900 400		61 7 3808 4871
E-mail	: tps@tps.com.au	Web:	www.tps.com.au